

### X-Cite exacte



# Stability. Control. Repeatability.

## Setting a new standard.

The X-Cite *exacte* has the convenience of pre-aligned, long-life lamps and easy installation combined with innovative functions in the newest generation of mercury arc fluorescence illuminators.

#### Stability You Can Count On

Using a DC-powered lamp coupled with proprietary Closed-Loop Feedback (CLF) technology, the X-Cite exacte maintains constant power and ensures ultra-stable illumination. The X-Cite exacte is ideal for long term time-lapse and comparative imaging, remaining consistent whether experiments last for milliseconds or days.

## Calibrated Output for Ultimate Repeatability

The X-Cite *exacte* can be calibrated using the X-Cite Optical Power Meter, an optional accessory. Ensure truly repeatable experiments by taking the variability out of illumination intensity.

#### Ideal for Live Cell Imaging

Equipped with an iris to provide adjustment of light intensity in steps as small as 1% from 0-100% the X-Cite *exacte* optimizes specimen illumination. This provides the flexibility and control of intensity attenuation required where photobleaching is a concern.

#### **High Speed Internal Shutter**

The X-Cite *exacte* is equipped with a high-speed internal shutter activated by push-button, foot pedal, RS-232 commands or TTL, eliminating the need to purchase additional hardware for high-speed applications.

The rapid response time of less than 6ms\* for time critical exposures protects delicate specimens.





A custom graphical user-interface (GUI) enables computer control of all X-Cite exacte functions, logs events and records and displays real-time stability data. Ideal for verifying intensity stability during automated or long term imaging experiments.



Designed with Safety in Mind

The X-Cite exacte has safety designed into all aspects of its operation: Interlock switches and a light guide detection system prevent accidental exposure to UV and intense visible light. Intelli-Lamp® monitors the system temperature to prevent overheating.

#### The X-Cite Advantage


All of the convenient and cost-saving features that made the original X-Cite a leader in fluorescence illumination are also designed into the X-Cite exacte. The pre-aligned Intelli-Lamp means easy set-up, and a 2000 hour lamp life guarantee means low operating costs. The X-Cite lamp provides superior excitation for a wide range of fluorophores across the entire spectrum. A variety of specially engineered adaptors ensures compatibility with microscopes from all major manufacturers.

The X-Cite exacte gives you convenience and control never before available. Engineered specifically to let you focus your energy where it belongs - on the best possible results.

For superior convenience and control, insist on X-Cite. Our work relies on accurate quantitative measurements of fluorescence intensities in cells over the duration of an experiment which can be a few hours or multiple weeks. The exacte provides both the short-term stability with the Closed-Loop Feedback and long-term stability with the calibrated light output. The exacte has greatly enhanced our capability to do quantitative fluorescence imaging work.

Damir Sudar, Staff Scientist Lawrence Berkeley Laboratory, Berkeley CA


#### **Spectral Comparison**

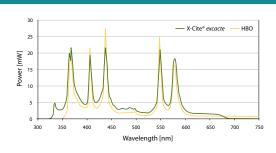

#### The Benefit of Closed-Loop Feedback

X-Cite exacte uses a Closed-Loop Feedback signal to monitor and stabilize intensity. As the lamp ages, the iris position is automatically adjusted to maintain intensity at the original set point.



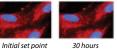





DC Lamp without Closed-Loop Feedback:








Initial set point

30 hours

60 hours 90 hours

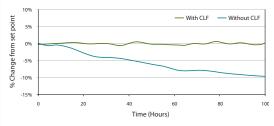

FEATURES	BENEFITS
DC lamp with Closed-Loop Feedback	Short and long-term intensity stability
Calibration via optional Optical Power Meter	Provides output in watts for consistent illumination and empirical repeatability
Adjustable Iris with 1% steps	Intensity control limits photobleaching
Internal High Speed Shutter	Flexibility in configuring imaging systems for different applications
Liquid Light Guide	Excellent illumination uniformity, no heat transferred to microscope
Automatic Light Guide Detection	Ensures proper/complete insertion of light guide to optimize performance and lifetime
Pre-aligned, 2000 hour lamp	Easy lamp replacement with no alignment required, lower running costs
Intelli-Lamp Technology	Maintains optimum lamp temperature, tracks lamp hours
Computer Control & GUI	Enables computer-driven applications and stability data exporting

TECHNICAL SPECIFICATIONS	
Includes	Lamp module, liquid light guide with adaptor, software CD, foot pedal, USB cable, grounded and shielded power cord, user guide, quick-start guide
Lamp	200W Mercury Vapor Short Arc, DC
Lamp Life	2000 hours guaranteed; 2500 hours typical
Adaptors	Compatible with microscopes from all major manufacturers
Filter	Bandpass, 340-675nm
Power Supply	High efficiency, switch mode, line isolated 100-120VAC / 200-240VAC, 50/60Hz
Warm-up Period	4-minutes (typical)
Connectivity	RS-232 control via USB port; PC and Mac control of lamp, iris shutter; TTL control of shutter via BNC; ready for integration in automated system
Drivers	Base commands same as X-Cite 120PC; Build onto existing driver scripts to control X-Cite exacte.
Panel Controls	Power on/off, Display Mode Selection, Intensity up/down adjustment, Closed-Loop Feedback on/off, Shutter open/close
Panel Display	Accumulated lamp hours, intensity setting, status messages
Panel Status Indicators	Lamp on, shutter open, display mode, calibration, light guide insertion, Closed-Loop Feedback
Dimensions (LxWxH)	13.5" x 7" x 8" 34cm x 18cm x 20cm
Weight	10.8lbs / 4.9kg
Certifications	CE marked, certified to IEC, Canadian and US Standards; RoHS compliant
Warranty	1 year (excluding lamp and light guide)
Patents	X-Cite exacte incorporates technology protected by the following patents: US# 5,521,392; US# 6,333,602; US# 6,437,861; US# 6,847,170; US# 7,335,901

Hg-LAMP CONTAINS MERCURY, Manage in Accord with Disposal Laws, see www.lamprecycle.org or +1 800 668-8752



www.excelitas.com x-cite@excelitas.com

L5N 6H7 CANADA Fax: +1 905 821-2055

2260 Argentia Road Telephone: +1 905 821-2600 Mississauga, Ontario Toll Free (USA and CAN): +1 800 668-8752